

# Detection and Cross Contamination Evaluation of Roundup Ready Soy via Mechanical Shear Homogenization using the GLH-850

Shari Garrett, Omni International, Inc.

## Introduction

Genetically modified (GM) crops are food products that have had their DNA modified to introduce new traits such as resistance to pests, diseases or environmental conditions. One popular GM crop is Roundup Ready® soy by Monsanto. Roundup Ready® soy is genetically engineered soybeans designed to be tolerant of the herbicide Roundup®. The Roundup® herbicide is a glyphosate that inhibits the enzyme 5-enolpyruvylshikimate-3-phosphate synthase (EPSPS) (1). EPSPS catalyzes an essential step in the biosynthesis of the amino acids phenylalanine, tyrosine and tryptophan in plants(2). When sprayed Roundup® quickly inhibits the growth of plants by interfering with the synthesis of these essential amino acids. Roundup® is typically used to inhibit the growth of weeds but usually at the cost of reduced crop yields. Round-up Ready® soybeans have been genetically engineered to express a version of the EPSPS gene that confers resistance to glyphosphates therefore allowing it to survive the doses required to kill weeds that infect farmers fields (1).

In order to determine the detection of Roundup Ready® soy, the soy products are homogenized and the DNA is purified from a variety of extraction techniques. Following extraction, the genetically modified EPSPS gene is detected via polymerase chain reaction (PCR). Rotor stator homogenizers are one of the best methods for processing a variety of samples from soybeans to soy powder. The GLH-850 is Omni's newest programmable rotor stator homogenizer. The GLH-850 can quickly process samples at speeds of

up to 25,000 RPM and sample volumes up to 10 liters. Particularly, the GLH-850 can process samples using a variety of stainless steel and Omni Tip plastic generator probes. Omni's disposable plastic generator probes can process a range of sample matrices and can eliminate cross contamination, samples loss and carryover concerns when dealing with sensitive PCR assays.

Herein, we evaluate the GLH-850's capability of processing a variety of soy sample matrices for the detection of Roundup Ready® soy. In conjunction, the potential for sample cross contamination from the use of a standard stainless steel generator probes was evaluated.

## Materials & Methods

### Equipment

- **General Laboratory Homogenizer (GLH) 850.** (CAT #20-010)
- **10 mm Stainless Steel Wide Window Generator** (Cat #G10-95W)
- **7 mm Hard Tissue Plastic Omni Tip** (Cat#30750H)
- **Omega Bio-Tek E.Z.N.Z Plant DNA Kit** (Cat#D3495-01)

7 mm Hard Tissue Plastic Omni Tip (Cat#30750H)



10 mm Stainless Steel Wide Window Generator (Cat #G10-95W)



## Sample Preparation

0.5 g of organic soybeans and 0.5 g of soy powder was processed in 2.5 ml of buffer CBL (Omega) with the GLH-850 fitted with a 10 mm stainless steel wide window generator probe at 20,000 RPM for approximately 15-30 seconds until a complete homogenate was achieved. The generator probe was disassembled and rinsed under a stream of DD H<sub>2</sub>O between samples. An additional 0.5 g of organic soybeans and 0.5 g of soy powder was processed in 2.5 ml of Buffer CBL with the GLH-850 fitted with a 7 mm Omni hard tissue plastic tip. The samples were processed at 20,000 RPM for up to 60 seconds until a complete homogenate was achieved. Each sample was processed with its own individual plastic generator probe.

### General Laboratory Homogenizer (GLH) 850.

(CAT #20-010)



## DNA Isolation, Quantification and Separation

500 µl of each homogenate was added

to 10 µl of 2-mercaptoethanol, vortexed and incubated at 65°C for up to 30 minutes (15 minutes for soybean). After incubation, the DNA was purified per the kit manufacturer's instructions. 1 µl of each eluant was analyzed for DNA quantification on the NanoDrop Spectrophotometer (Thermo Fisher Scientific).

The size and quality of the isolated DNA was determined by agarose gel electrophoresis. Approximately 300 ng of DNA was mixed with 10 µl of TBE/Urea loading dye (BioRad) and separated on a 1% agarose gel in TBE running buffer (BioRad). Electrophoresis was carried out at 140V for 45 minutes. The gel was stained in ethidium bromide for 15 minutes and visualized on a GelDoc EZ System (images/data not show).

### PCR Assay and Fragment Analysis

Primers for the amplification of the 5'-enolpyruvylshikimate-3-phosphate synthase (EPSPS) gene were chosen using published sequences. Primers were synthesized by Integrated DNA Technologies (Iowa, USA).

**Table 1. PCR Primers**

Name	Sequence (5'-3')	Reference
RRO1	TGGCGCCCAAAGCTTGCATGCGC	3
RRO4	CCCCAAGTTCCTAAATCTTCAAGT	3

Each amplification reaction was established by adding 1x reaction buffer; 10 mM Tris-HCl, 50 mM KCl, 1.5 mM MgCl<sub>2</sub>, 0.2 mM dNTPs, 5% Glycerol, 25 U Taq (New England Biolabs), 0.5 µM of each primer and 20 ng of template. Amplification was carried out using the T100 Thermal Cycler by BioRad per the settings in table 2 (4).

**Table 2. PCR Program**

	Temperature	Time
Hot Start/Denaturation	95°C	3 minutes
50 Cycles	95°C	45 seconds
	60°C	45 seconds
	72°C	25 seconds
Final Extension	72°C	10 minutes

PCR products were analyzed on a 2% agarose gel and stained using ethidium bromide. Fragments were visualized and analyzed via BioRad's Gel Doc EZ system.

**Table 3. Soy Bean and Soy Powder DNA concentrations**

Sample	Stainless Steel Generator Probe	Omni Plastic Generator Probe
Soybean	174.9 ng/µl	114.0 ng/µl
Soy Powder	284.9 ng/µl	217.0 ng/µl

### Results


In this study, we investigated the GLH-850's capability of homogenizing a variety of soy matrices for the detection of Roundup Ready® resistance genes. In conjunction, we observed the potential for cross contamination when using stainless steel generator probes. When performing sensitive assays such as PCR, cross contamination can lead to erroneous and false positive results. In this particular assay, soybeans and soy powder was processed using the GLH-850 using both stainless steel and plastic generator probes. Genomic DNA was extracted from each sample and the glyphosate resistance EPSPS gene was detected via end-point PCR.

#### Genomic DNA Quantification

The isolated genomic DNA was quantified by spectrophotometry. DNA yields averaged 144.5 ng/µl and 250.9 ng/µl of soybean and soy powder respectively (Table 3).

#### PCR Detection


The fragment length of each amplicon from each sample type was analyzed on a 2% agarose gel. The expected size of the amplified EPSPS gene is shown to be 356 bp. As shown, the soy powder samples contained bands at the expected base pair length for the EPSPS gene (Figure 1).

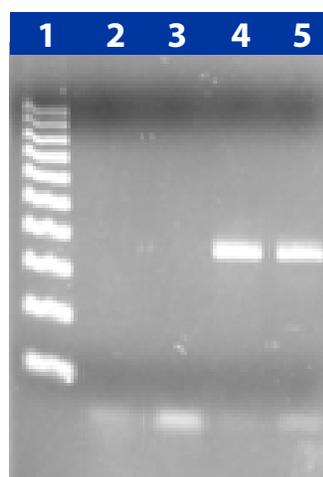

**Figure 1** Roundup Ready Detection from Disposable Omni Tip. Lane 1: 100 bp Ladder, 2-3: Soybean, 4-5: soy powder

The Roundup Ready® resistance gene was not detected for the organic soybean samples while a single fragment at the desired length was detected from the soy powder samples.

However, when processing with a single stainless steel generator probe cross contamination was evident in the soybean sample. Despite disassembling and rinsing the probe, trace amounts of soy powder DNA remained on the probe and tainted the soybean sample creating a false positive (Figure 2). In order to eliminate the chance of cross contamination when using a stainless steel generator probe it is imperative that the probe is disassembled, thoroughly cleaned and autoclaved.

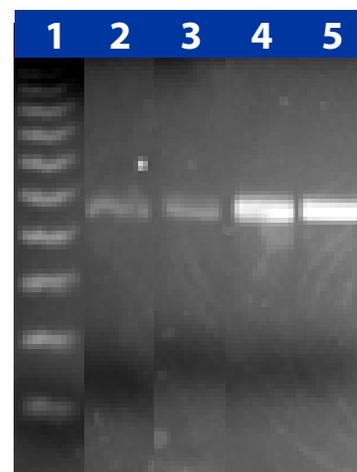

**Figure 2** Roundup Ready Detection from Stainless Steel Probe. Lane 1: 100 bp Ladder, 2-3: Soybean, 4-5: Soy Powder

### Conclusion

Rotor stator homogenizers such as the GLH 850 are capable of disrupting a variety of sample matrices for DNA extraction with subsequent PCR analysis. The GLH 850 is excellent at processing soy samples for the extraction of genomic DNA in excess of 110 ng/µl. Using either the stainless steel or plastic generator probes the DNA was shown to be of excellent quality for PCR analysis. However, it is noteworthy that when processing multiple samples with a stainless steel generator probe carryover is very likely without the use of

thorough cleaning techniques. In cases where very sensitive assays are performed, Omni's plastic Omni Tip generator probes are highly recommend since they can be disposed of after each sample. This is particularly important when GM foods are being detected and testing needs to accurate and reliable.

Roundup® and Roundup Ready® are registered trademarks of Monsanto Technology LLC

## References

1. Funke, T., Han, H., Healy-Fried, M. L., Fischer, M., & Schonbrunn, E. (2006). Molecular basis for the herbicide resistance of Roundup Ready crops. *Proceedings of the National Academy of Sciences*, 103(35), 13010-13015.  
Retrieved February 29, 2016.
2. Tzin, V., & Galili, G. (2010). The Biosynthetic Pathways for Shikimate and Aromatic Amino Acids in *Arabidopsis thaliana*. *The Arabidopsis Book / American Society of Plant Biologists*, 8, e0132. <http://doi.org/10.1199/tab.0132>
3. Studer, E., Rhyner, C., Luthy, J., & Hubner, P. (1998). Quantitative competitive PCR for the detection of genetically modified soybean and maize. *Zeitschrift Für Lebensmitteluntersuchung Und -Forschung A*, 207(3), 207-213.  
Retrieved February 29, 2016.
4. Tengel, C. (2001). PCR-Base Detection of Genetically Modified Soybean and Maize in Raw and Highly Processed Foodstuffs. *BioTechniques*, 31(2), 426-429.  
Retrieved February 29, 2016.



935-C Cobb Place Blvd. NW  
Kennesaw, GA 30144  
800.776.4431 • 770.421.0058  
[www.omni-inc.com](http://www.omni-inc.com)

